

- 1. CHARGE: VOLTAGE DIAL @ 0
 TURN AC ON FIRST FOR 2 min.
 TO ALLOW VACCUM TUBES TO
 WARM UP. THEN DC SWITCH ON.
- 2. THEN TURN VOLTAGE CONTROL TO NO MORE THAN 175. WAIT 3 min. TO CHANGE (AMPS TO 0). THEN TURN DC OFF.
- 3. DISCHARGE WITH ROD.
 DISCHARGE AGAIN.
 DISCHARGE AGAIN.
 THEN RETURN TO @ REST.